

***The H105 ultrasonic cleaner has an internal clock.
Here is how to set it...***

The screen is touch sensitive and options can be chosen in one of two ways:

- Press and hold your selection until selected
- Press and quickly release then press and release **Select** at the bottom of the screen

From the home screen select **Settings**

From the Main menu select **Settings**

From the Settings menu select **Settings 2**

From the Settings 2 menu select **Time & date**

To select and change the Year, Month, Date or time, move up and down the menu by pressing **Up** and **Down** or press on **'Year'**, **'Month'** or **'Day'** to select.

When Year, Month or Day is highlighted pressing **Select** will increase the value. Press and hold to cycle through the values.

Once the year, month and day are correct you may set the time.

Select Time from the menu. Press **-** or **+** to decrease or increase the time.

Press **Set** to save. Return to the home screen by pressing **Back** until it is displayed.

If you have any problems either email sales@walkerelectronics.co.uk or contact a member of our team on 01636 892410.